3M Food Safety

3M[™] Petrifilm[™] Plates & 3M[™] Petrifilm[™] Plate Reader



Technician Productivity Maximized



3M[™] Petrifilm[™] Plates... Take a Closer




Count Plate

- Enumeration of Aerobic Total Flora.
- Reading facilitated by an indicator which colours colonies red.




Petrifilm Enterobacteriaceae Count Plate

- Enumeration of Enterobacteriaceae.
- Production of acid by colonies, highlighted through a pH indicator which turns yellow.
- Upper film traps the gas produced by glucose fermentation.

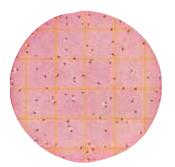

Petrifilm Coliform Count Plate

- Enumeration of coliforms (total or thermotolerant).
- Reading facilitated by an indicator which colours colonies red.
- Upper film traps the gas produced by lactose fermentation.

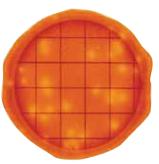

Petrifilm Rapid Coliform **Count Plate**

- Rapid enumeration of coliforms.
- Very sensitive pH indicator reveals production of acid by growing colonies.
- Growth acceleration reduces the response time.




Petrifilm E. coli /Coliform **Count Plate**

- Two tests in one: enumeration of Escherichia
- coli and coliforms. Contains a ß-glucuronidase indicator for detection of E. coli.
 - Results in 24 to 48 hours.



Petrifilm High-Sensitivity **Coliform Count Plate**

- Enumeration of small numbers of coliforms (less than 1 or 2 per gram).Inoculation of 5ml of sample
- or dilution of it.
- Reading advantages (colour and gas) identical to those for Petrifilm Coliform count plates.

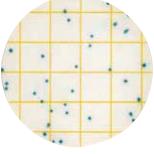

Petrifilm Select E. coli **Count Plate**


- Specific test for E. coli, only *E. coli* colonies are visible. Contains a ß-glucurcnidase
- indicator.






Petrifilm Yeast and Mold Count Plate

- Enumeration of yeast and mould population. Requires no addition of
- antibiotics.

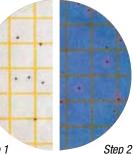

Step 1 Petrifilm Staph Express **Count System**

- Enumeration of
- Staphylococcus aureus. One or two step analysis
- (Dnase reaction for the
- second step).Results in 24 to 27 hours.

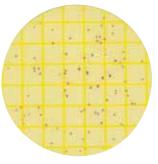


Petrifilm Environmental Listeria Plate

- Detection of Listeria in envi-. ronmental samples.
- Provides Quantitative, Semi-Quantitative and Qualitative results in 26-30 hours.
- No enrichment required.




Put 3M[™] Petrifilm[™] Plate technology to work for you.

3M Petrifilm[™] Count Plates: Coating technology

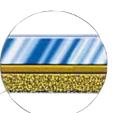

1. Top Film

Plastic Film coated with adhesive, indicator and cold water soluble gel.

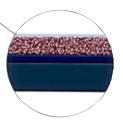

2. Bottom Film

Plastic coated paper printed with a grid, adhesive standard methods nutrients, cold water soluble gel.

Protecting companies and brands

While maximizing productivity is essential in today's business world, providing the highest quality product is paramount. You can trust Petrifilm Plates to reduce variability, from technician to technician, shift to shift, plant to plant.

More than 200 evaluations have been completed from validating agencies (AFNOR, NORDVAL, AOAC International, AOAC RI, EMMAS) and peer-review publications worldwide on the precise, consistent results delivered by Petrifilm Plates.' These are just some of the reasons Petrifilm Plates are in use by over 3000 companies in Europe.'


Three Simple Steps to improve analysis with 3M[™] Petrifilm[™] Plates



. Inoculate Petrifilm" Plates are easily inoculated. No media preparation is required.



 Incubate Petrifilm[®] Plates can be incubated in a smaller incubator.

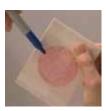

3. Interpret Simply count colonies. Indicator dyes facilitate interpretation.

Data on file.

3M[™] Petrifilm[™] Plates Benefits



Better working conditions for technicians

- Sample ready for immediate handling
- Eliminate media preparation time, routine tests done in half the time.
- Reduce time pressure, improve productivity
- Reduce burden on day-to-day laboratory logistics
- Allow more flexibility to organize laboratory activity
- Reduce control requirements for accreditated laboratory (a quality certificate is provided with each box)
- Standardized manufacturing process delivers consistent plates.

Compact Film Size

- Save space in the laboratory, improve working conditions
- Less storage capacity required, better inventory management

The Quality Standard from 3M

- Guaranteed manufacture from 3M
- Manufacturing site certified to ISO 9002 standard
- Strict quality control procedure helps reduce media variation

3M[™] Petrifilm[™] Plate Reader... **CONSISTENTLY ACCURATE COUNTS IN JUST 4 SECONDS.**

Automated plate reading is fast and simple

This compact machine fits efficiently into the lab environment and saves critical time by automating your lab's labor intensive plate reading. Technicians are free to tackle the other jobs that demand their attention, confident that the Petrifilm Plate Reader will handle its tasks with excellent performance, time after time. It processes the five most common plate tests (Aerobic Count, Coliform Count, E. coli/Coliform Count, E. coli Count and Enterobacteriaceae Count) to handle the majority of your plate reading needs. The Petrifilm Plate Reader:

- Reduces the chance for human error
- Decreases variations (enhances repeatability and reproduceability)
- Allows data management for audit purposes or FDA 21CFR Part11 compliance
- Can store actual color images of plates for further interpretation



Results and a color image of the plate appear on your computer screen, simplifying colony verification.



Log file data is automatically stored and easily transferred into a text file, making exporting of essential information easy.





3M[™] Petrifilm[™] Plate Reader

| 3M ID | Description | Packaging | Item No. | Plate Qty |
|----------------|------------------------------|-----------|----------|-----------|
| FH-5000-0164-6 | Petrifilm Plate Reader | 1 | 6499 | 1 |
| 78-8131-2876-2 | Petrifilm Plate Reader Stand | 1 | 6499S | 1 |

3M[™] Petrifilm[™]

| 3M ID | Description | Packaging | Item No. | Plate Qty |
|-------------|--|-----------|----------|-----------|
| 70200572157 | Detrifilm Aerobie count Diete | 20 x 50 | 6406 | 1000 |
| 70200572124 | Petrifilm Aerobic count Plate | 2 x 50 | 6400 | 100 |
| 70200572165 | Petrifilm Coliform count Plate | 40 x 25 | 6416 | 1000 |
| 70200572140 | | 2 x 25 | 6410 | 50 |
| 70200590142 | Petrifilm E. coli count Plate | 20 x 25 | 6414 | 500 |
| 70200572132 | | 2 x 25 | 6404 | 50 |
| 70200718859 | | 20 x 25 | 6435 | 500 |
| 70200718859 | Petrifilm Select E. coli count Plate | 2 x 25 | 6434 | 50 |
| 70200590159 | Detrifilm Veget (Maylel gount Diete | 20 x 50 | 6417 | 1000 |
| 70200561242 | Petrifilm Yeast/Mould count Plate | 2 x 50 | 6407 | 100 |
| 70200703885 | Detrifilm Llich Constituity Coliform Disto | 20 x 25 | 6415 | 500 |
| 70200703877 | Petrifilm High-Sensitivity Coliform Plate | 2 x 25 | 6405 | 50 |
| 70200647538 | Petrifilm High-Sensitivity Coliform Spreader | 2 | 6481 | 2 |
| 70200670936 | Petrifilm Enterobacteriacae count Plate | 40 x 25 | 6421 | 1000 |
| 70200670928 | Petrimini Emeropacteriacae count Plate | 2 x 25 | 6420 | 50 |
| 70200659525 | Detrifilm Danid Caliform count Dista | 20 x 25 | 6412 | 500 |
| 70200659535 | Petrifilm Rapid Coliform count Plate | 2 x 25 | 6402 | 50 |
| 70200734534 | Detrifilm Stoph Everage count Plate | 2 x 25 | 6490 | 50 |
| 70200734542 | Petrifilm Staph Express count Plate | 20 x 25 | 6491 | 500 |
| 70200734559 | Petrifilm Staph Express Disk | 20 | 6492 | 20 |
| 70200734567 | reumini Stapii Express Disk | 5 x 20 | 6493 | 100 |
| 70200719063 | Petrifilm Staph Express Spreader | 2 | 6425 | 2 |
| 70200742990 | Detrifilm Environmental Listoria Plata | 2 x 25 | 6447 | 50 |
| 70200743006 | Petrifilm Environmental Listeria Plate | 8 x 25 | 6448 | 200 |
| 70200743022 | Petrifilm Environmental Listeria Spreader | 2 | 6498 | 2 |



3M Health Care 3M Food Safety Europe, Middel East & Africa Carl-Schurz-Str. 1 41453 Neuss, Germany Phone: +49 (0) 2131/14-3000 Fax: +49 (0) 2131/14-4397

www.3M.com/microbiology